


Thank you for choosing INSPECTIS®

This product has been tested with Windows 7/10/11 Professional. Below you will find step by step instructions on how to install.

If your device requires Windows drives, install the drivers before running **INSPECTIS®**. You can find device drivers on the provided memory stick together with other installation files.

Following Inspectis devices require driver installations:

Item Nr.	Name
HD-130	Inspectis 4K UHD PCIe Capture Device
DM-008 / IN-001	Inspectis Digital Microscope, Capillaroscope

1. Install the software by running "**INSPECTIS x.x Setup**" installation file. Follow on-screen instructions.
2. Once installation is finished, right click on INSPECTIS icon  on your computer desktop and select **Run as Administrator** first time.
2. Start Inspectis. On "Select Inspectis Version" dialog box, select:
 - a. "**INSPECTIS® ProX**" if you have ordered a **ProX** license. Insert the provided software protection USB dongle into a free USB port of your computer. Enter license number and other requested information. Follow up on-screen instructions to complete installation.
 - b. "**INSPECTIS® Pro**" if you have ordered a **Pro** license. Insert the provided software protection USB dongle into a free USB port of your computer. Enter license number and other requested information. Follow up on-screen instructions to complete installation.
 - c. "**INSPECTIS® Basics**" if you have ordered a **Basics** license. Enter license number and other requested information. Follow up on-screen instructions to complete installation.
 - d. "**INSPECTIS® OAI**" if you have ordered a **OAI** license. Enter license number and other requested information. Follow up on-screen instructions to complete installation.
 - e. 15days trial. All software functions will be enabled for 15 days. License number is not required.
3. Software will be registered and activated automatically over internet once you provide license number and other requested information. If on-line registrations is prohibited or not possible due to your network restrictions, complete the Off-line Registration form and email it to info@inspect-is-com to get an activation key.
4. Select your video capture adapter or any other Inspectis compatible video device and software will start monitoring live image of the digital microscope.

For more information refer to instruction's manual of **INSPECTIS®** software.